

# ELISA Assay for Glutathione S-Transferase Alpha

For Research Use Only

# INTRODUCTION

Glutathione S-Transferase (GST) has multiple isoforms. This assay is specific for Glutathione S-Transferase Alpha (GSTA) and is not known to cross react with the mu, pi, or theta variants. GSTA is a common biomarker for hepatocellular damage.<sup>1</sup> It also conjugates GSH to 4-hydroxynonenal, a product of lipid peroxidation<sup>2</sup> and is an important player in cellular antioxidant defense mechanisms.<sup>3</sup>

## **PRINCIPLES OF PROCEDURE**

This is a standard sandwich enzyme-linked immunosorbent assay (ELISA). The plate is pre-coated with anti-GSTA and blocked, ready for the addition of samples and standards. The assay should take approximately 3 hours to run, plus any required sample preparation time.

# MATERIALS PROVIDED

Component	Description	Volume	Storage	Cat no.
Anti-GSTA Plate	96-well microplate coated and blocked	1 plate	4°C	GS41a
Assay Buffer	Buffer used to dilute samples and reagents	100 mL	4°C	GS41b
10x Wash Buffer	Buffer used to wash the plate	30 mL	4°C	GS41c
GST Alpha Standard	10 μg/mL GSTA	20 µL	4°C	GS41d
<b>Detection Antibody</b>	Anti-Human-GSTA	130 µL	4°C	GS41e
HRP-Conjugate	Streptavidin-HRP conjugate	130 µL	4°C	GS41f
TMB Substrate	Stabilized TMB color reagent	20 mL	4°C	GS41g

# MATERIALS NEEDED BUT NOT PROVIDED

- 1. Microplate reader with 450 nm filter
- 2. Adjustable micropipettes and tips
- 3. 3N Sulfuric Acid (H<sub>2</sub>SO<sub>4</sub>)
- 4. Deionized Water (dH<sub>2</sub>O)

## STORAGE

- 1. Store the components of this kit at the temperatures specified on the labels.
- 2. Unopened reagents are stable until the indicated kit expiration date.

# WARNINGS AND PRECAUTIONS

- 1. Use aseptic technique when opening and dispensing reagents.
- 2. This kit is designed to work properly as provided and instructed. Additions, deletions or substitutions to the procedure or reagents are not recommended, as they may be detrimental to the assay.

GST Alpha ELISA Product Number: GS41 Store at 4°C FOR RESEARCH USE ONLY Control Number: GS41.210817 Page 1 of 4

#### **PROCEDURAL NOTES**

- 1. Reagents can be used immediately upon removal from refrigeration.
- 2. To minimize errors in absorbance measurements due to handling, wipe the exterior bottom of the microplate wells with a lint-free paper towel prior to inserting into the plate reader.
- 3. Do not save excess or diluted reagents for future use.

## SAMPLE STORAGE

Samples should be stored at  $-80^{\circ}$ C and thawed just prior to use. Avoid repeated freeze/thaw cycles for best results. This assay was developed and validated with human serum samples, however it does cross-react with rat.

#### SAMPLE PREPARATION

It is recommended to do multiple sample dilutions to ensure that the concentration falls within the accepted range for the assay. Samples should be assayed neat or diluted 1:2 in Assay Buffer.

## **REAGENT PREPARATION**

- 1. **GST Alpha Standard:** Immediately prior to use, dilute 1:1000 by adding 10 μL of Standard to 10 mL of Assay Buffer, giving a final concentration of 10 ng/mL.
- 2. **10x Wash Buffer:** Dilute the wash buffer 1:10 by adding 30 mL of 10x Wash Buffer to 270 mL of dH<sub>2</sub>O.
- 3. **Detection Antibody:** Immediately prior to use, dilute 1:100 by adding 120 µL of Detection Antibody to 12 mL of Assay Buffer.
- 4. **HRP-Conjugate:** Immediately prior to use, dilute 1:200 by adding 60 μL of HRP-Conjugate to 12 mL of Assay Buffer.

#### **STANDARD CURVE PREPARATION**

Set up for the standard curve by labeling dilution tubes and dispensing the indicated volumes of Assay Buffer and 10 ng/mL Standard Stock Solution according to Table 1 below.

Standard	GST Concentration (ng/mL)	Assay Buffer (µL)	Volume of 10 ng/mL Standard (µL)	Final Volume (µL)		
S7	10	-	1000	1000		
S6	8.0	200	800	1000		
S5	5.0	500	500	1000		
S4	2.5	750	250	1000		
S3	1.0	900	100	1000		
S2	0.5	950	50	1000		
<b>S</b> <sub>1</sub>	0.25	975	25	1000		
S <sub>0</sub>	0	1000	-	1000		

**Table 1:** Standard Curve Preparation

## ASSAY PROCEDURE

- 1. Add 100  $\mu$ L of Standards and Samples to the corresponding wells on the microplate in duplicate. Incubate at room temperature for one hour. See Scheme 1 below for a suggested plate layout.
- 2. Dump the contents of the plate and wash each well three times with 300 µL of Wash Buffer. After the final wash, tap the plate on a lint-free paper towel to make sure there is no solution left in the wells.
- 3. Add 100 µL of the Detection Antibody to each well. Incubate at room temperature for one hour.
- 4. Wash the plate as in step 2.
- 5. Add 100 µL of the HRP Conjugate to each well. Incubate at room temperature for 30 minutes.
- 6. Wash the plate as in step 2.
- 7. Add 100  $\mu$ L of TMB Substrate to each well. Allow the color to develop for 30 minutes at room temperature.
- 8. Stop the reaction by adding 25  $\mu$ L per well of 3N Sulfuric Acid (H<sub>2</sub>SO<sub>4</sub>).
- 9. Read the plate at 450 nm in a microplate reader.


Scheme 1. Suggested I fate Layout												
	1	2	3	4	5	6	7	8	9	10	11	12
Α	s <sub>0</sub>	s <sub>0</sub>	U <sub>1</sub>	Ul	U9	U9	U17	U17	U25	U25	U33	U33
В	$s_1$	$s_1$	U <sub>2</sub>	U <sub>2</sub>	U10	U10	U18	U18	U26	U26	U34	U34
С	s <sub>2</sub>	s <sub>2</sub>	U3	U3	U11	U11	U19	U19	U27	U27	U35	U35
D	s <sub>3</sub>	s <sub>3</sub>	U4	U4	U12	U12	U20	U20	U28	U28	U36	U36
Е	S4	S4	U5	U5	U13	U13	U21	U21	U29	U29	U37	U37
F	S5	S5	U <sub>6</sub>	U <sub>6</sub>	U14	U14	U22	U22	U30	U30	U38	U38
G	S <sub>6</sub>	S6	U7	U7	U15	U15	U23	U23	U31	U31	U39	U39
Н	S7	S7	U8	U8	U16	U16	U24	U24	U32	U32	U40	U40

Scheme 1: Suggested Plate Layout

# CALCULATIONS

- 1. Average all duplicate well absorbance values.
- 2. Subtract the average absorbance values for the blank wells (S<sub>0</sub>) from all other well pairs.
- 3. Plot a standard curve using the corrected absorbance values of each Standard (y-axis) versus the Standard concentration (x-axis).
- 4. Determine the concentration of each unknown using the equation of the line.

## Figure 1: Typical Standard Curve

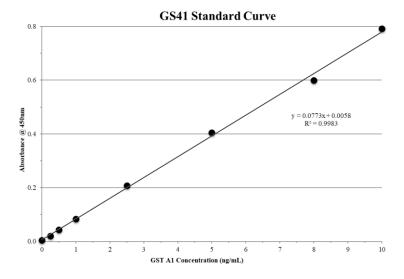

#### REFERENCES

- 1. Vaubourdolle, M, et al.; (1995) Clinical Chemistry 41:1716-1719
- 2. Awasthi, Y. C, et al.; (2004) Free Radic. Biol. Med. 37: 607-619
- 3. Yang, Y., et al.; (2002) Toxicol. Appl. Pharmacol. 182: 105-115

#### DISCLAIMER

This information is believed to be correct but does not purport to be all-inclusive and shall be used only as a guide. Oxford Biomedical Research, Inc. shall not be held liable for any damage resulting from handling or from contact with the above product. See catalog for additional terms and conditions of sale.

#### **ORDERING INFORMATION**

For additional kits or a complete catalog please visit our website at www.oxfordbiomed.com or call 800-692-4633.

#### **TECHNICAL SUPPORT**

If you need technical information or assistance with assay procedures, call our Technical Support Department at info@oxfordbiomed.com, or 800-692-4633. Our staff will be happy to answer your questions about this or any other product in the Oxford Biomedical line.

#### **GUARANTEE AND LIMITATION OF REMEDY**

Oxford Biomedical Research, Inc. makes no guarantee of any kind, expressed or implied, which extends beyond the description of the materials in this kit, except that these materials and this kit will meet our specifications at the time of delivery. Buyer's remedy and Oxford Biomedical Research, Inc.'s sole liability hereunder is limited to, at Oxford Biomedical Research, Inc.'s sole liability hereunder is limited to, at Oxford Biomedical Research, Inc.'s sole liability hereunder is limited to, at Oxford Biomedical Research, Inc.'s option, refund of the purchase price of, or the replacement of, material that does not meet our specification. By acceptance of our products, Buyer indemnifies and holds Oxford Biomedical Research, Inc. harmless against, assumes all liability for the consequence of its use or misuse by the Buyer, its employees, or others. Said refund or replacement is conditioned on Buyer notifying Oxford Biomedical Research, Inc. within thirty (30) days of the receipt of product. Failure of Buyer to give said notice within thirty (30) days of receipt of product shall constitute a waiver by the Buyer of all claims hereunder with respect to said material(s).

Oxford Biomedical Research, Inc. P.O. Box 522 Oxford, MI 48371 U.S.A.

Orders: 800-692-4633 Technical Service: 248-852-8815 Fax: 248-852-4466 E-mail: <u>info@oxfordbiomed.com</u>

Made in the U.S.A.